



NanoFlowSizer



This project has received funding from the European Union's Horizon 2020 research and innovation program under grant agreement No 855867.

A novel instrument with new capabilities for in-line & offline sizing of nanosuspensions

**Rut Besseling
(Scientific Director)**

on behalf of the InProcess-LSP team

www.inprocess-lsp.com






Process Analytical Technology (PAT) solutions




Webinar outline

1. Challenges in nanosuspension (Process) Analysis
2. Spatially Resolved DLS and the benefits
3. Application examples and considerations
4. Conclusions

1. Challenges in (Process) Analysis of nanosuspensions

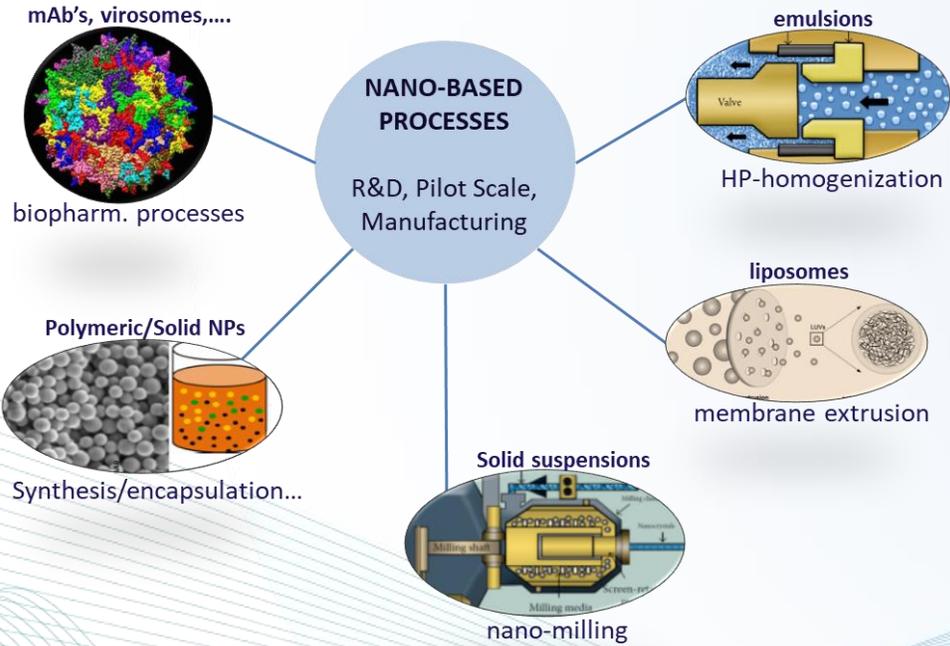



- Strong growth in R&D & Manufacturing of nanosuspensions
- Precise & high quality Manufacturing (Pharma/Chemical/....)
- Required Process/Product knowledge => Process Analytics

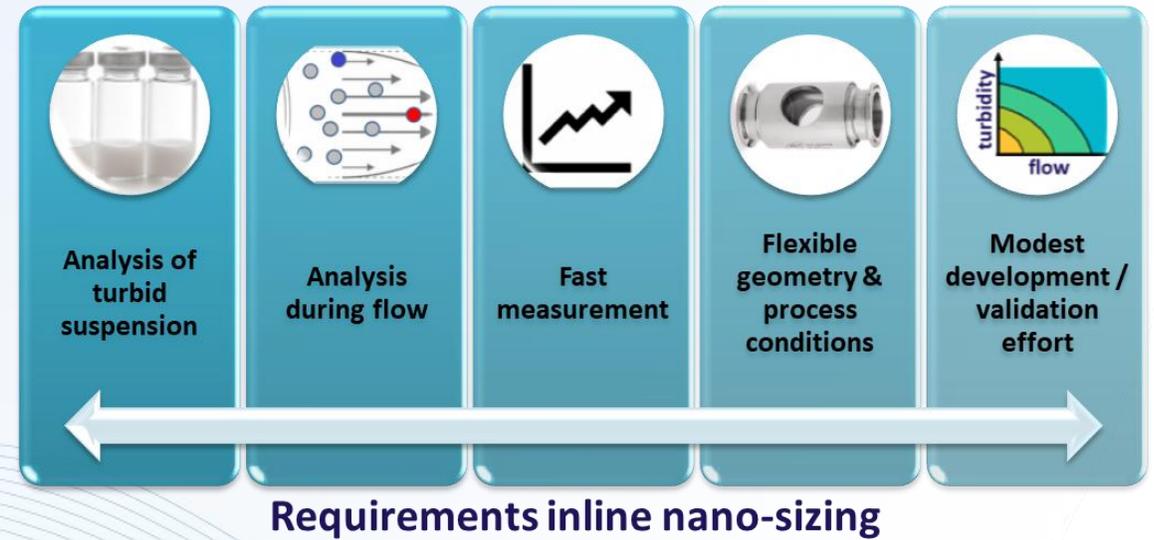





Noninvasive/PAT measurement needs beyond 'standard' instrument capabilities

1. New approach: Spatially Resolved DLS

- ❖ **'Low coherence' Dynamic Light Scattering** with 'instant' depth resolution:
 - Measurement of highly turbid samples by pathlength selection
 - Measurement in process/flow conditions:
 - resolve local flow & extract Brownian diffusion
- ❖ **Rapid process monitoring:** 10s per measurement
- ❖ **Flexible and fully non-invasive implementation possibilities**
- ❖ NIR-source (1300nm): optimal for highly turbid samples
- ❖ 180° backscattering: limits effect of large particles in polydisperse samples.

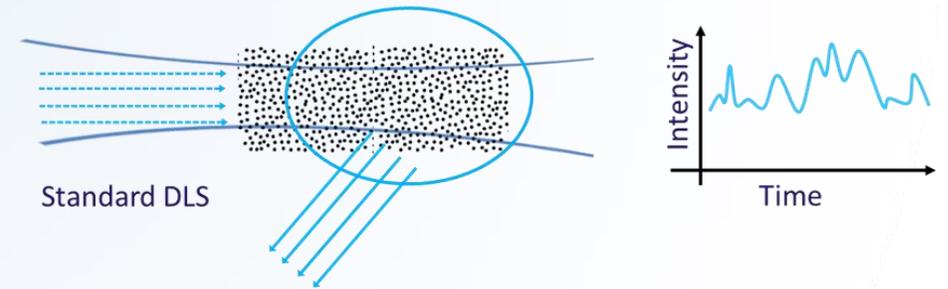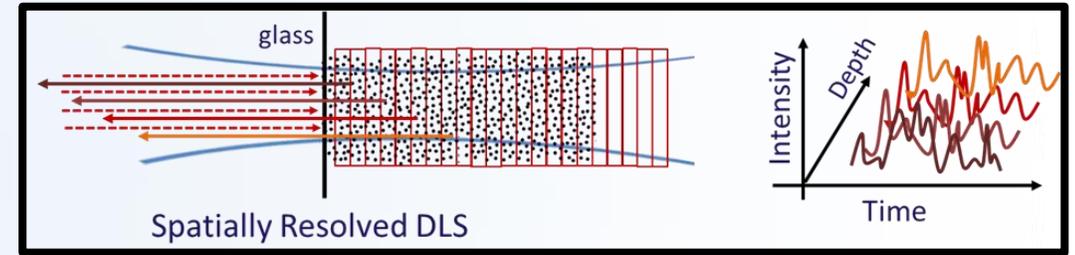


New opportunities for Industrial and Academic R&D

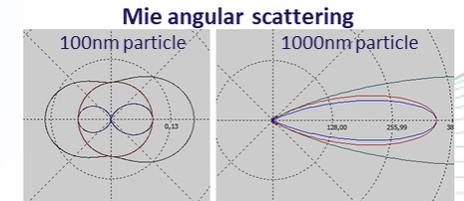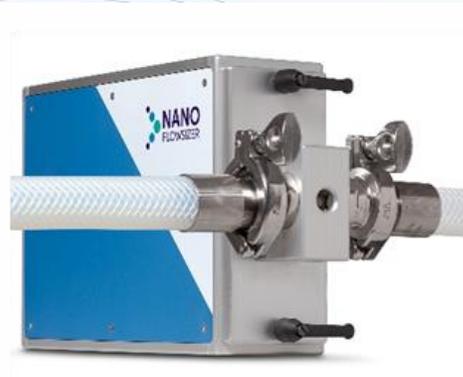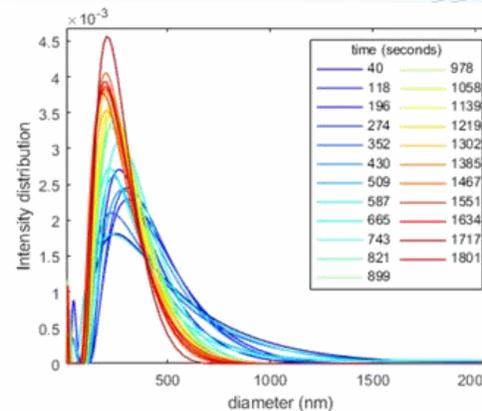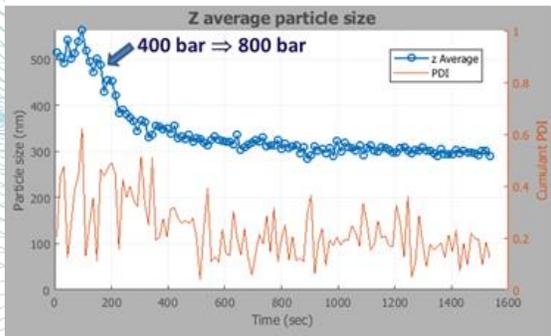

2. Background: standard Dynamic Light Scattering

standard DLS records fluctuations in scattered laser intensity due to Brownian diffusion


⇒ fluctuation rates via signal correlation ⇒ Diffusion constant(s) ⇒ hydrodynamic size (distribution)

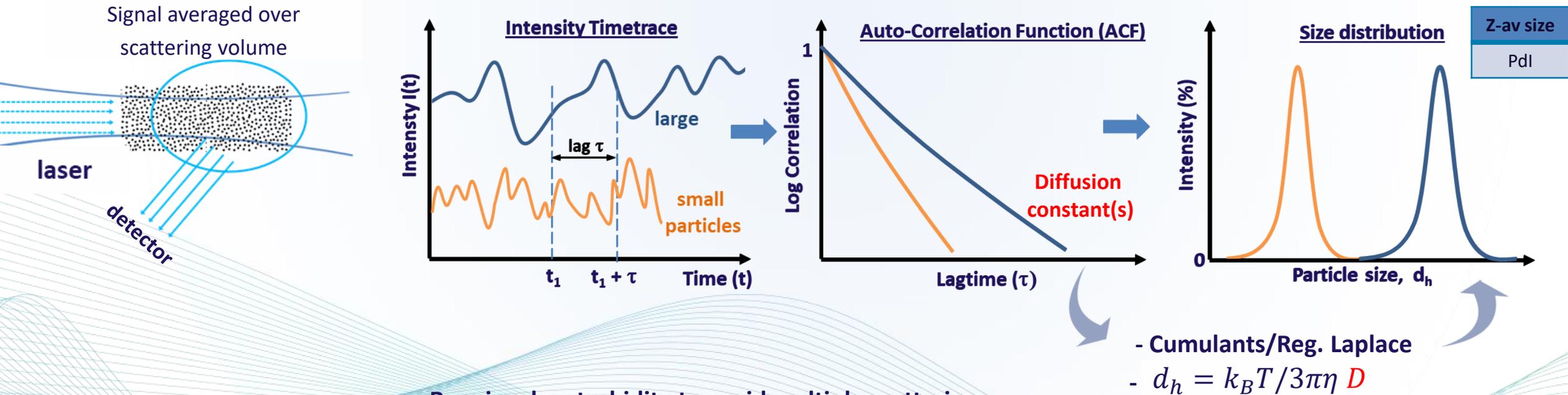


- Requires low turbidity to avoid multiple scattering
- Requires absence of flow

2. Spatially Resolved Dynamic Light Scattering


add the spatial dimension to standard DLS:

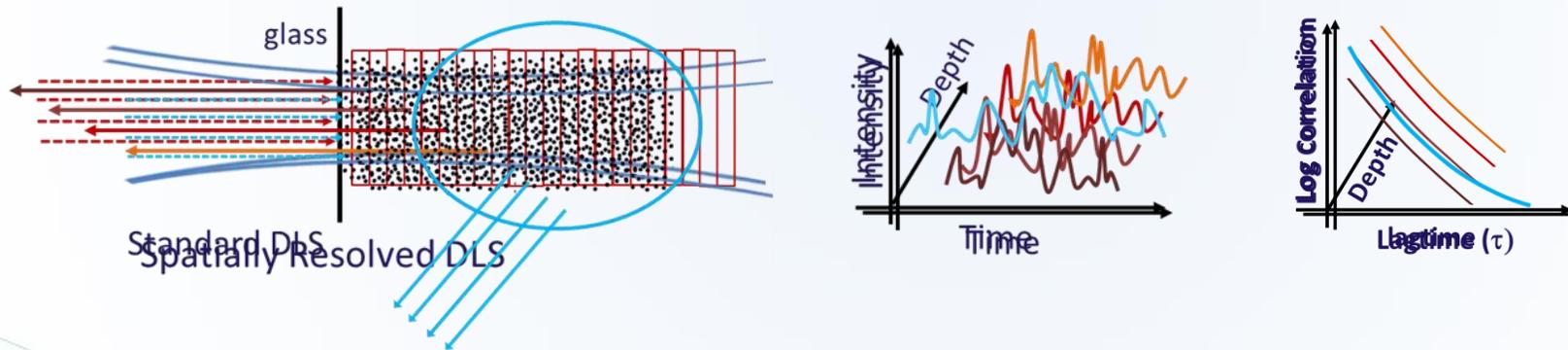

2. Spatial information via Low Coherence Interferometry

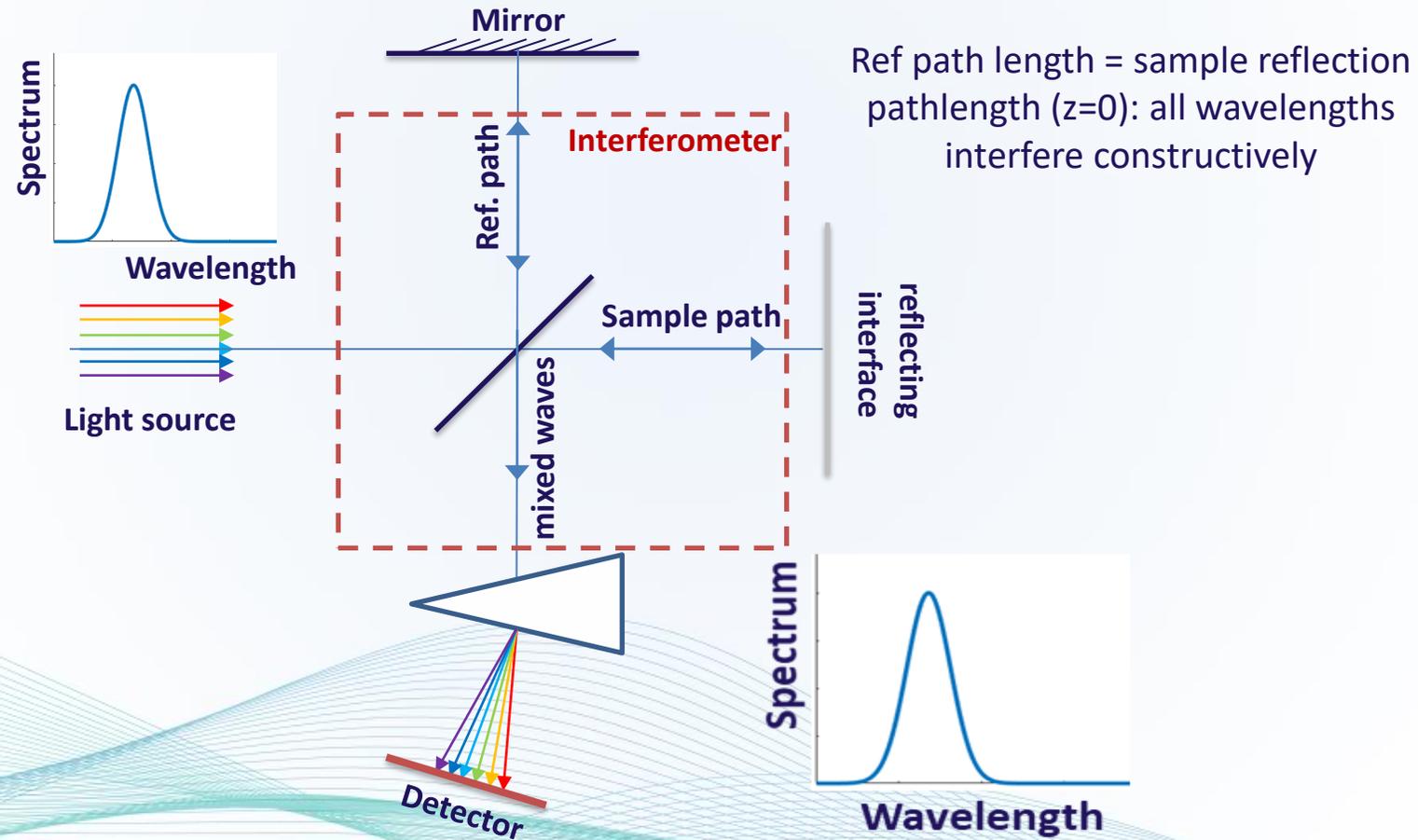


2. Spatial information via Low Coherence Interferometry

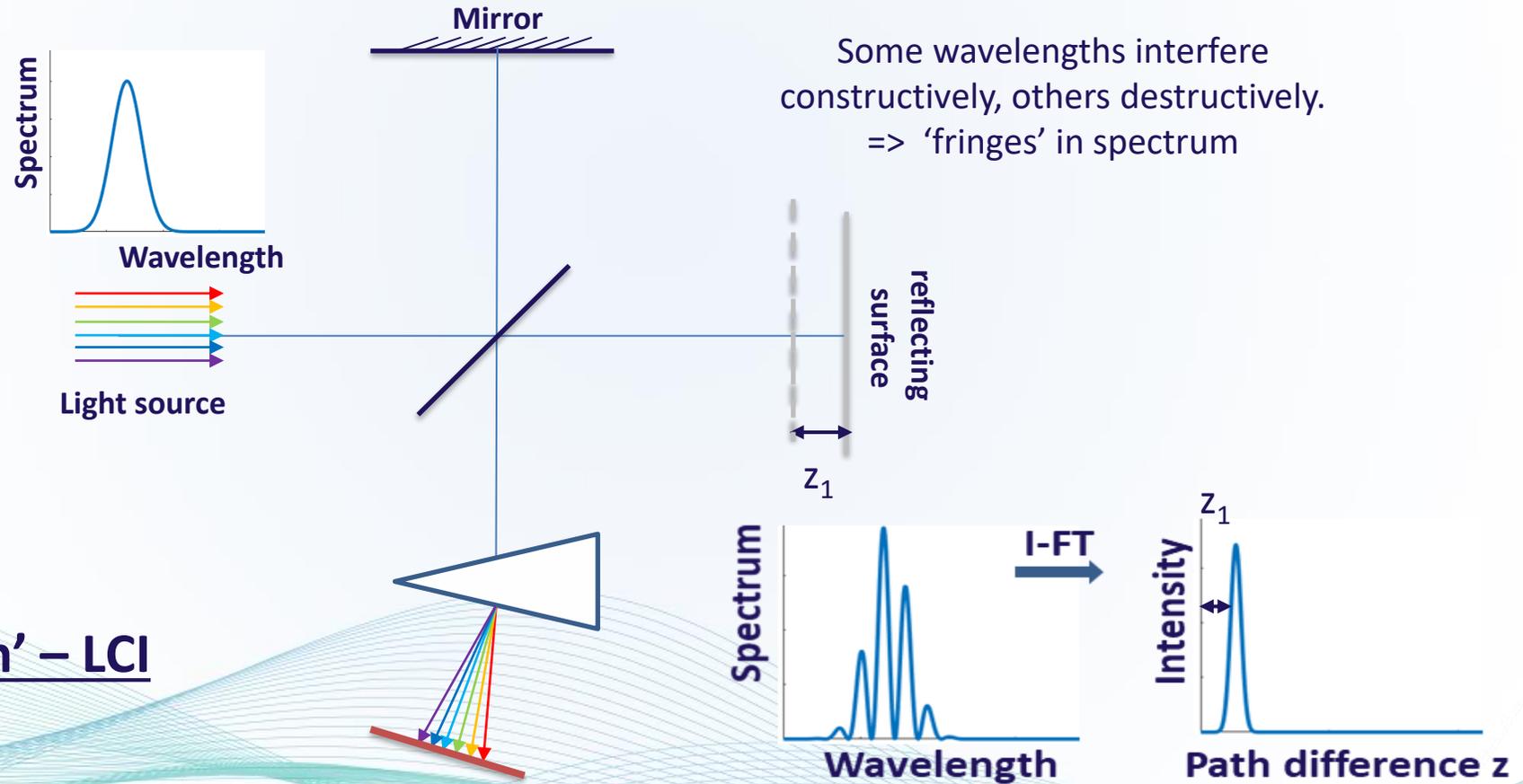


'Fourier-Domain' – LCI

2. Spatial information via Low Coherence Interferometry

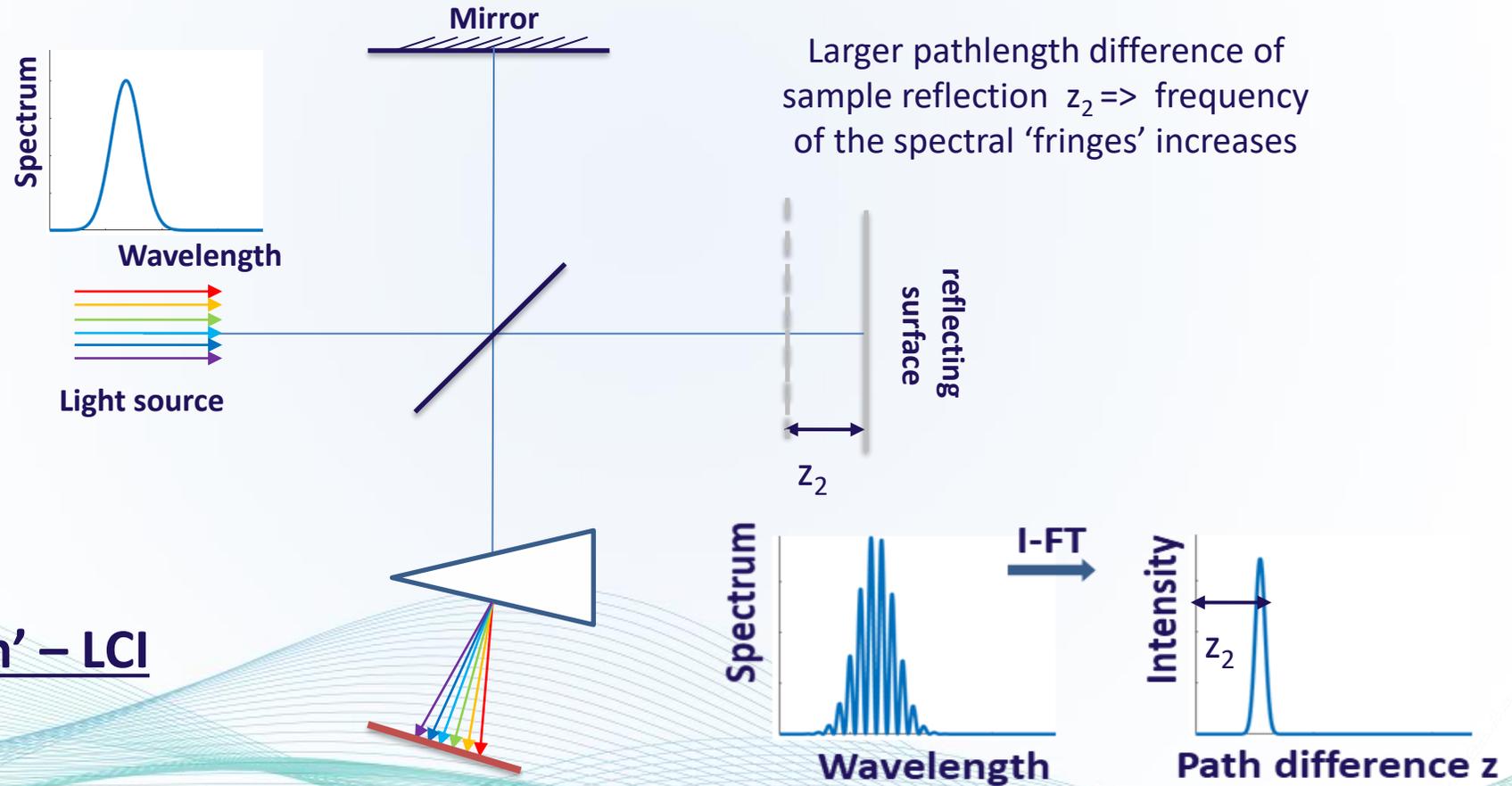


2. Spatial information via Low Coherence Interferometry

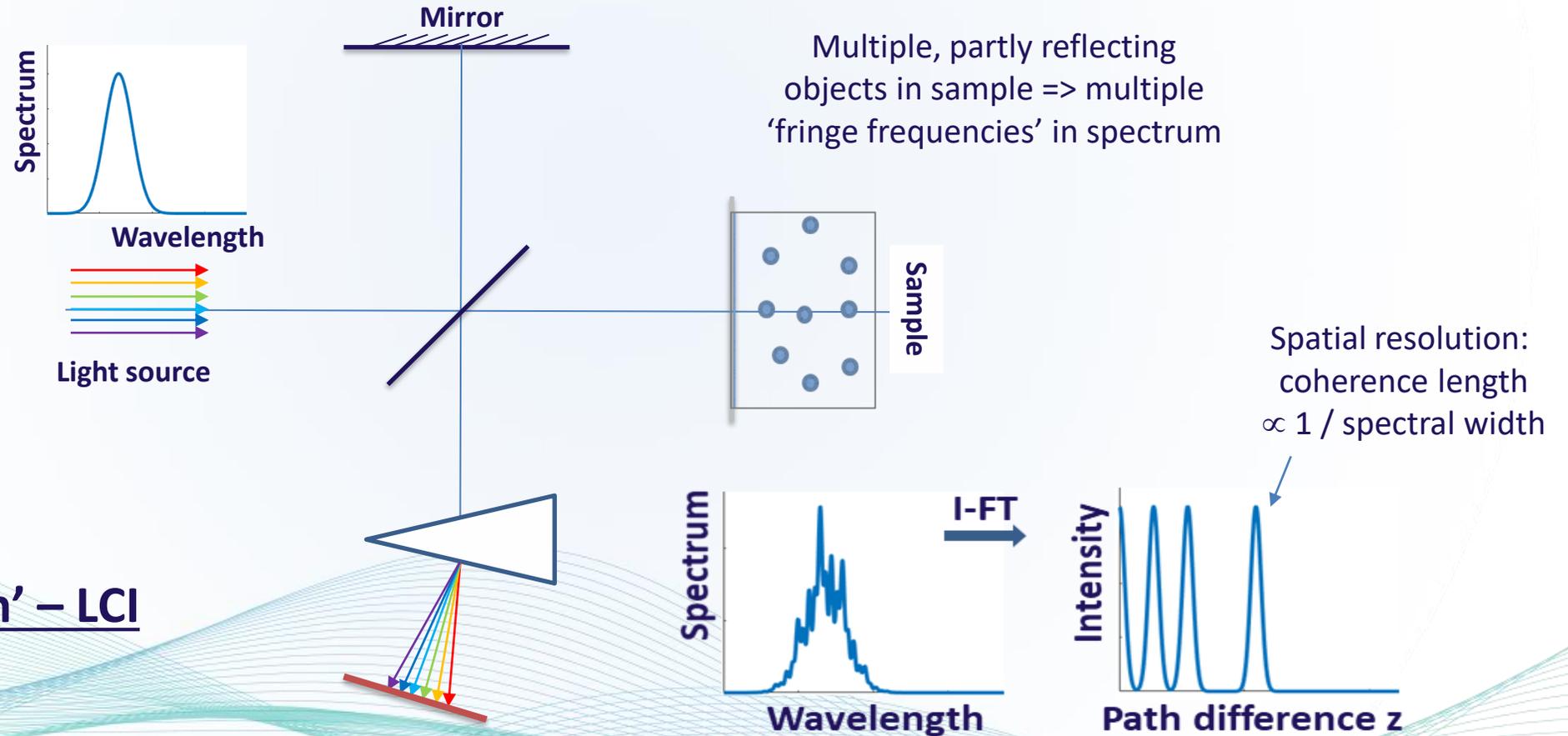




'Fourier-Domain' – LCI

2. SR-DLS in practice

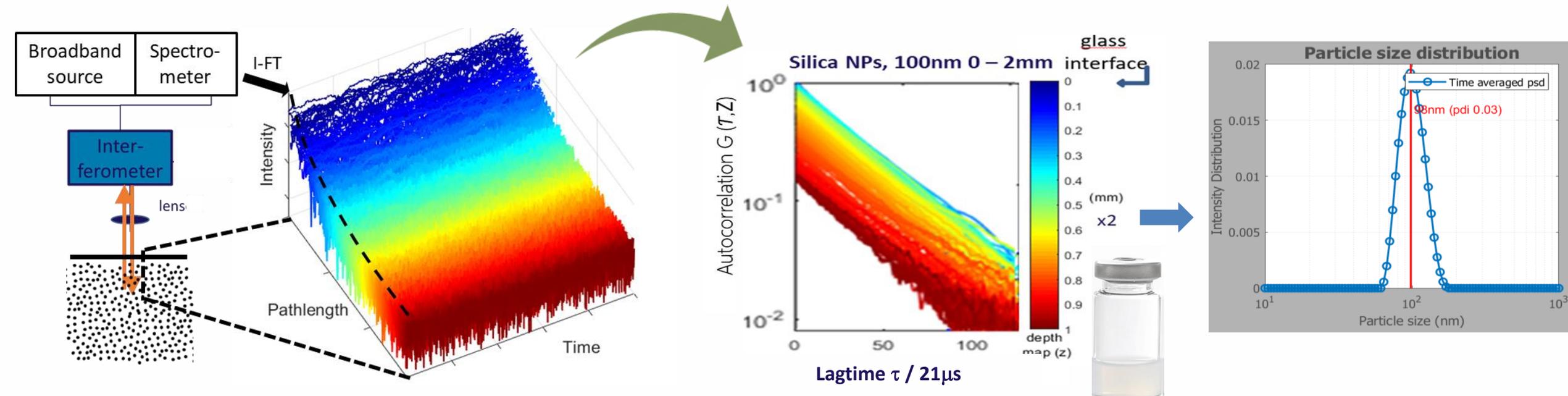

- Acquisition rate 76kHz \Rightarrow $I(z)$ resolved 'instantly' \Rightarrow Fluctuations $I(t, z)$ resolved at each depth.
- > 1000 ACF's per measurement \Rightarrow Resolve particle motions over $\sim 2.5\text{mm}$
- High information content \Rightarrow Z-av and Pdl in few seconds

2. SR-DLS benefits: high turbidity

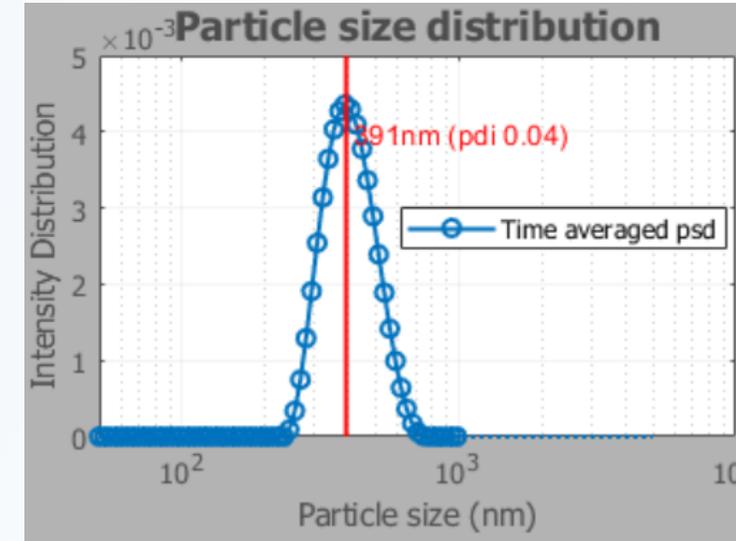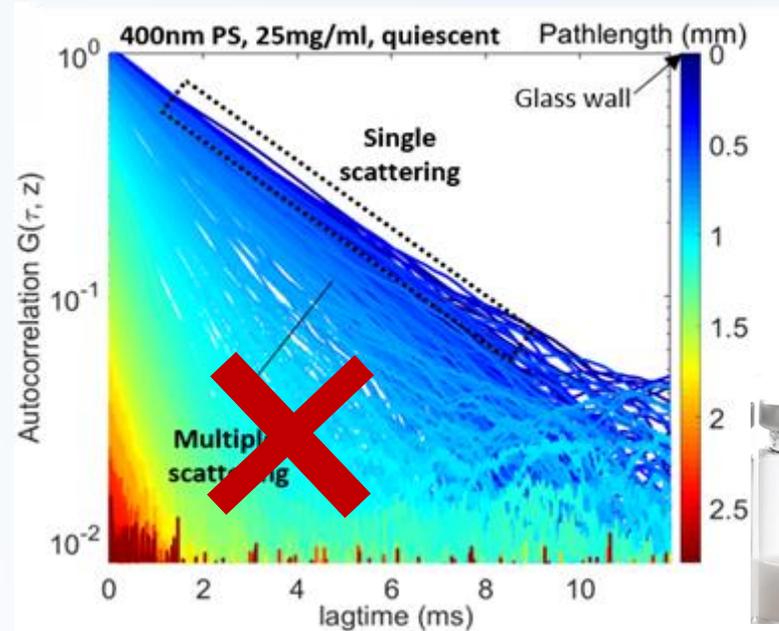
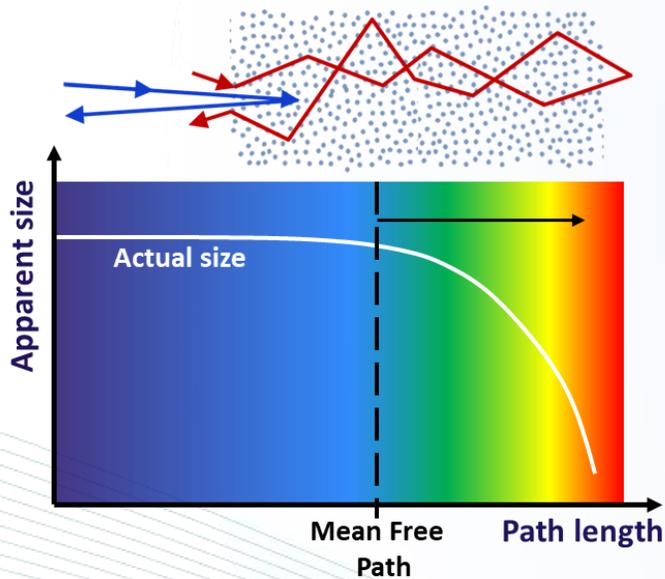


- Single & Multiple scattering regimes diagnosed automatically
- Multiple scattering data filtered (SR-DLS)
- Advantage: no dilution needed for turbid samples *


* photon mean free path down to $\sim 100\mu\text{m}$ @1300nm
e.g. Polystyrene suspensions up to at least 5 vol% (size dependent)

2. SR-DLS benefits: measurement in flow

Depth

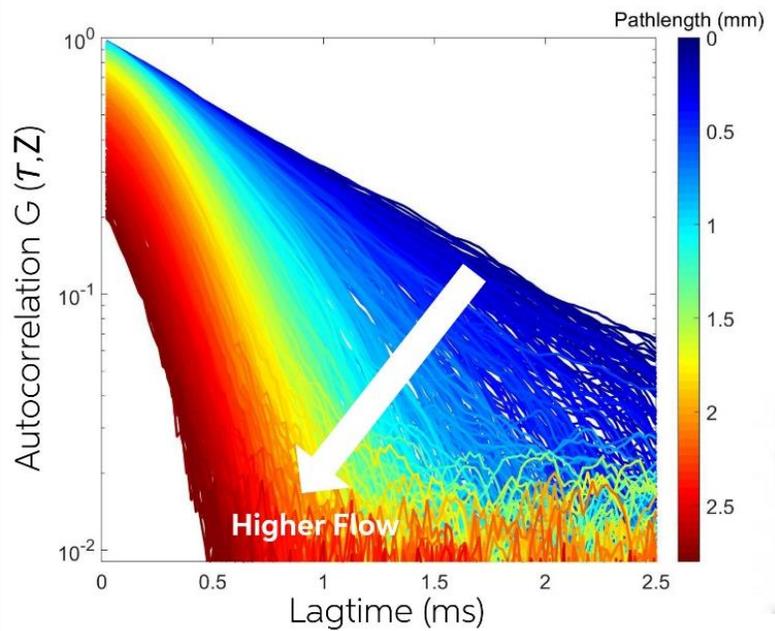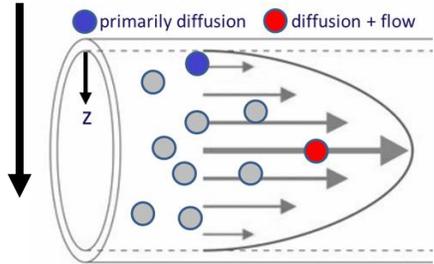

∅=22mm

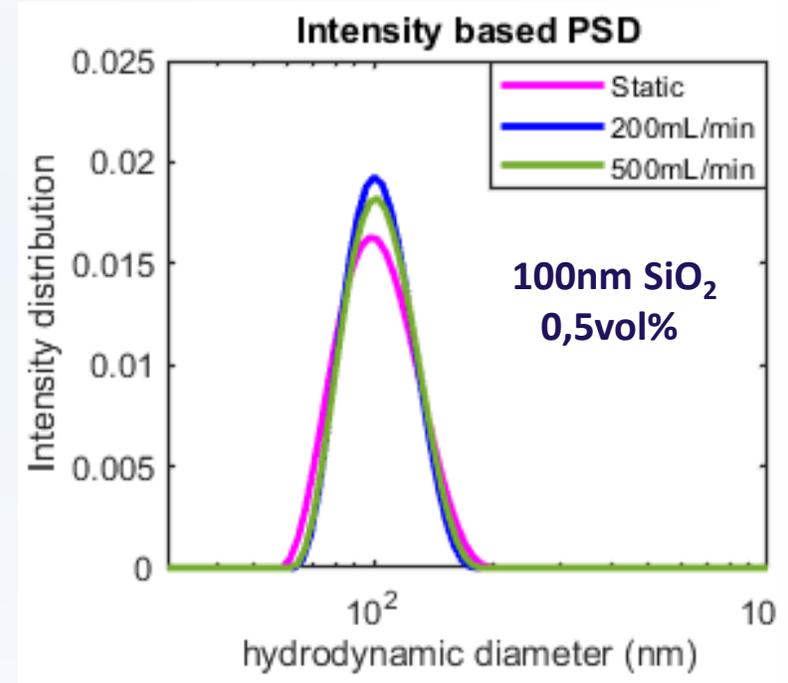

Diffusion




$$G(\tau, z) = G_{v=0}(\tau) \times G_{\text{flow}}(\tau/\tau_{\text{flow}}(z))$$

beam passage time

- ACF's contain information on diffusion and local flow speed.
- Flow measured and corrected for => size analysis from DIFFUSIVE PART OF THE ACF
- Flow-independent size measurement. No prior knowledge of flow required.

2. NanoFlowSizer system

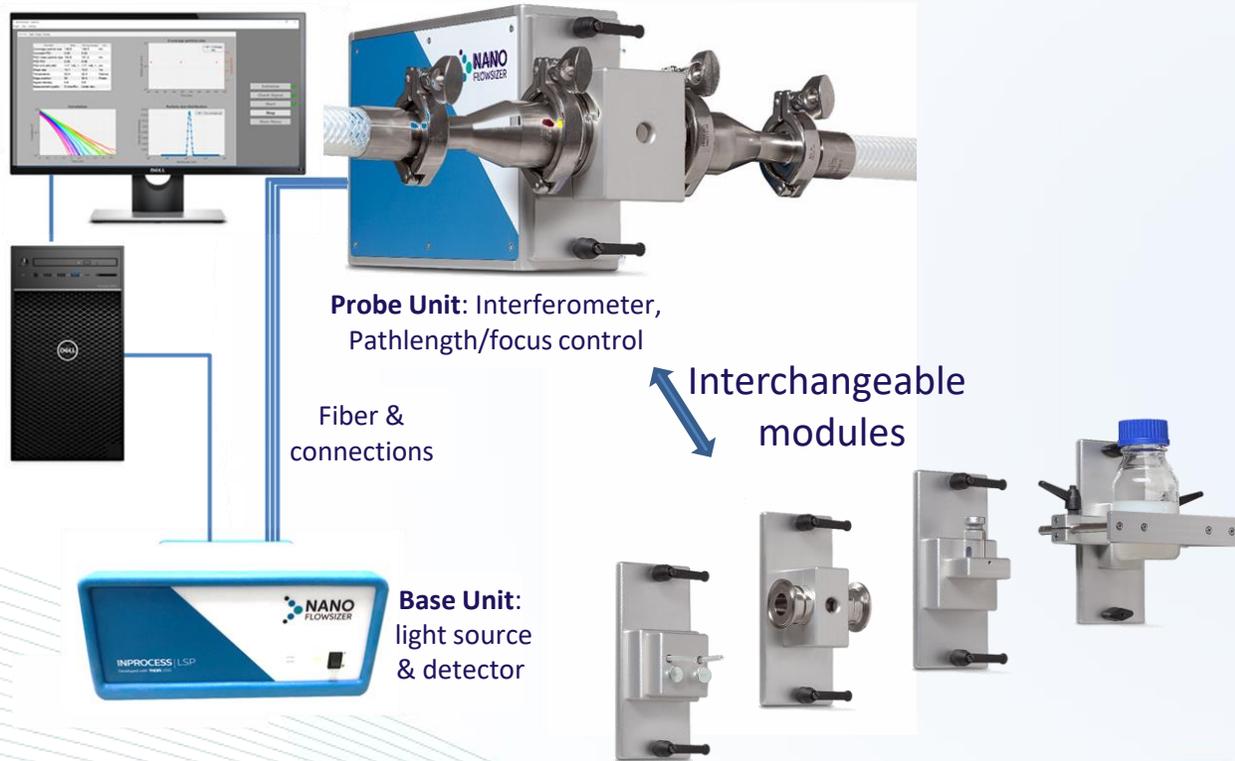

XsperGo software:




- handling ~Gb/s data in a continuous manner
- Real time measurement (~10s per point)

Outputs:



- Z-average, PDI (ISO22412)
- PSD: Intensity/volume, $d_{10,50,90}$
- Correlation function
- Flow and Scatter Profile
- Data Quality Diagnostics
- Advanced time & depth resolved diagnostics

MEASUREMENT SPECIFICATIONS	
Central wavelength	1300 nm
Axial scan rate	Up to 76 kHz
Maximum measuring depth / optical path length	3.5 mm
Maximum scattering volume	5 nL
Detection	Backscattering (180°)
Depth resolution	4 μm
Particle size range	10 - 1000 nm (at least)
Maximum process flow	6 L/min for 100nm polystyrene (5 mg/ml) in water


3. Applications: μ -flow cell & TiO₂ synthesis

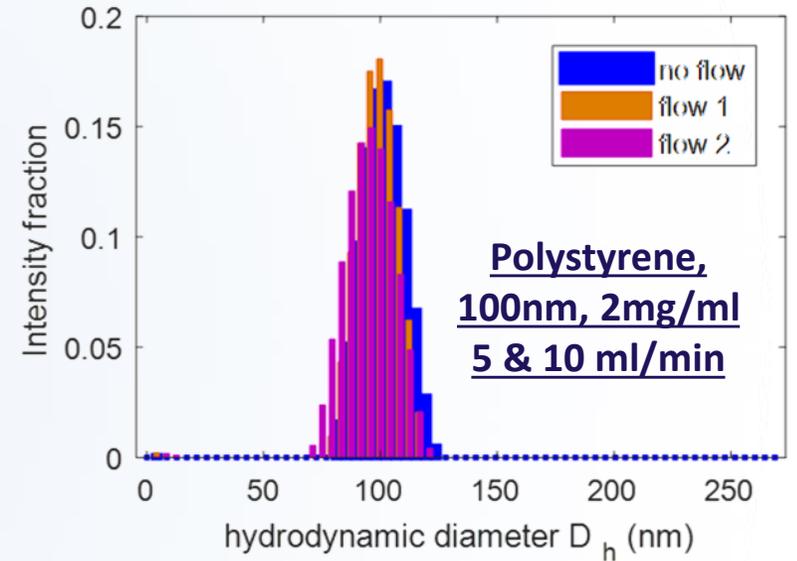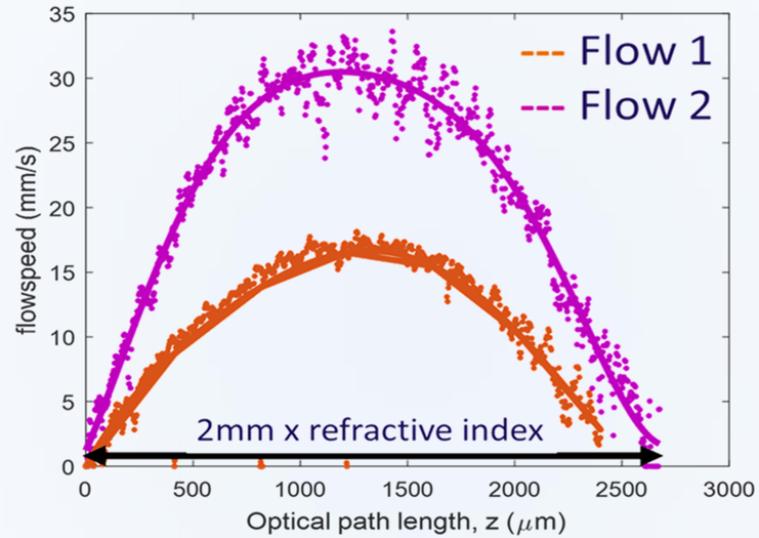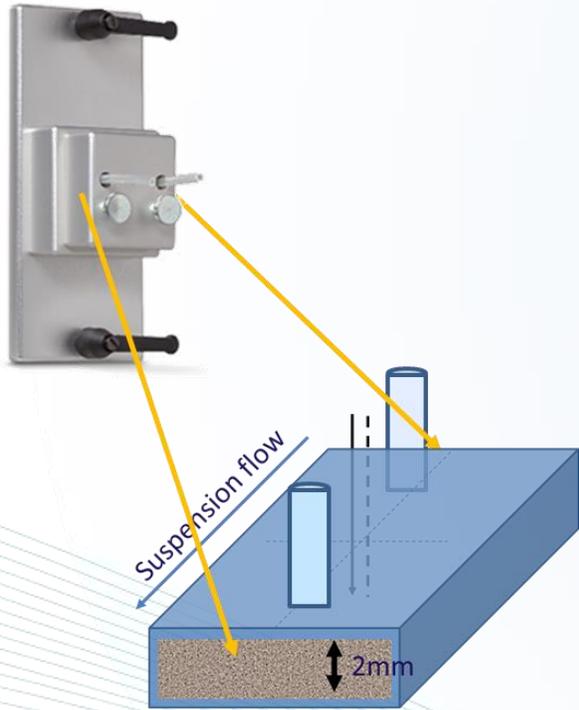



TiO₂ synthesis microflow sampling loop

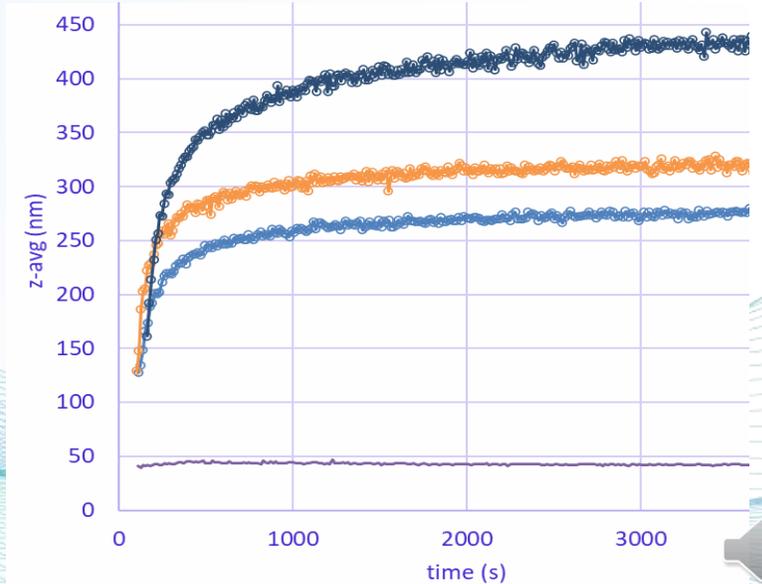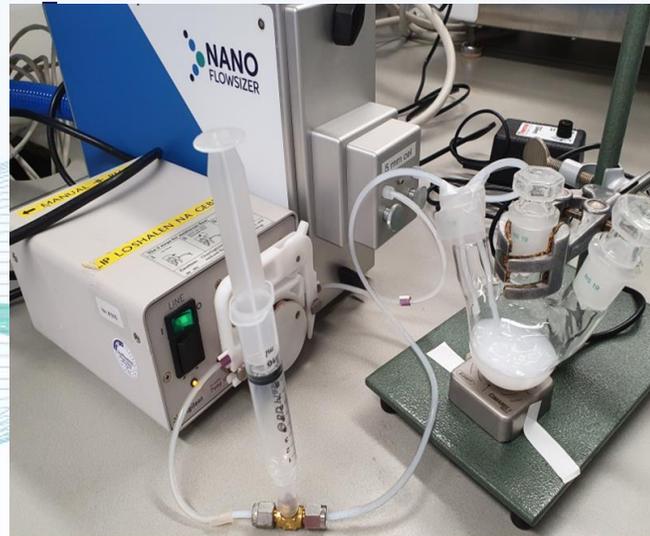

3. Monitoring SiO₂ synthesis & aggregation

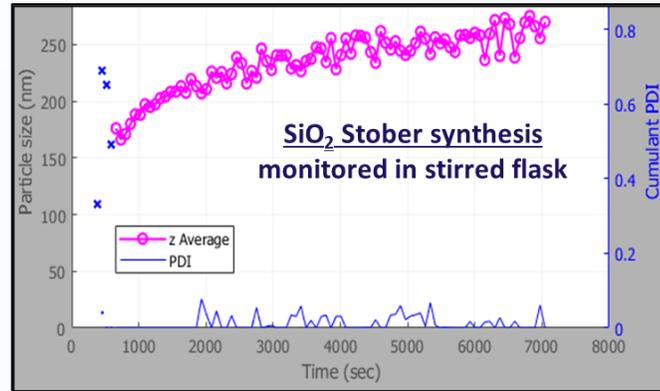

+ measured scatter intensity →




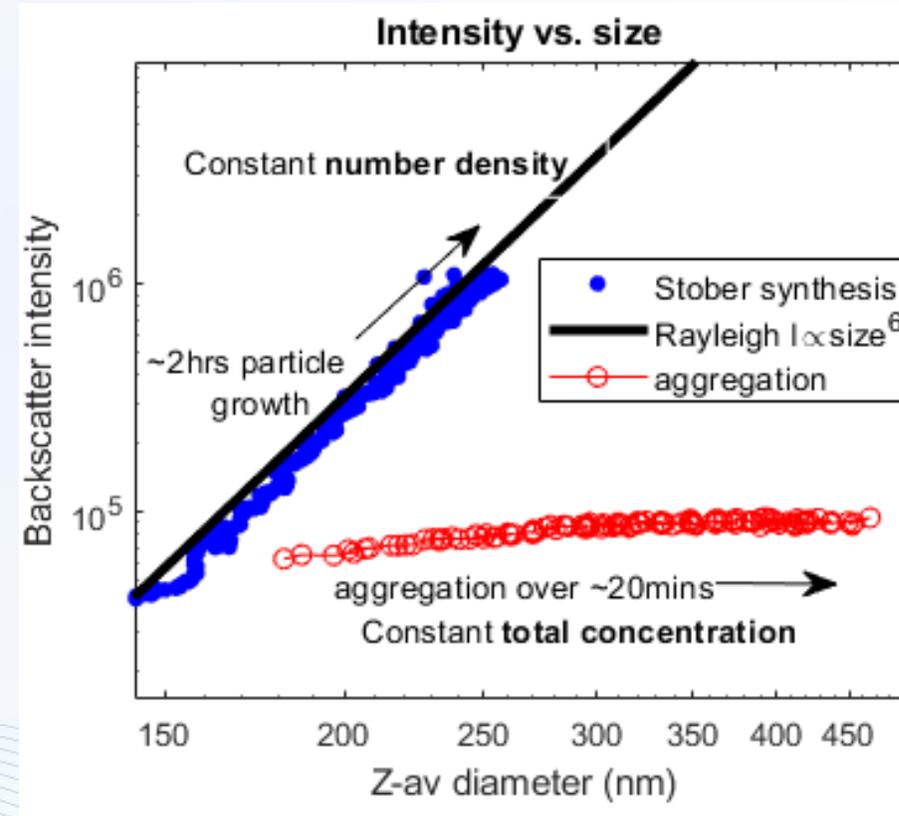
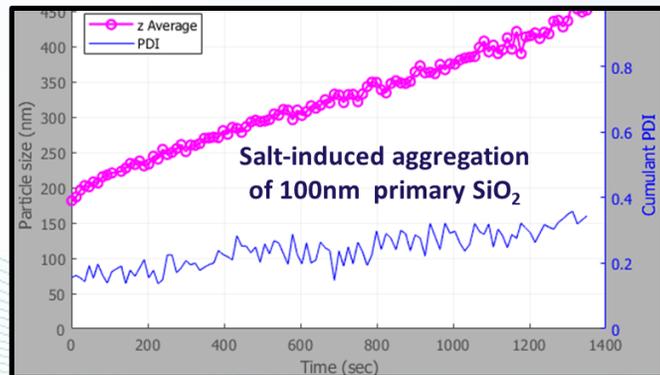


Note:
 $\lambda_{NFS} = 1300\text{nm}$
 Rayleigh law valid
 up to $d \sim 250\text{nm}$

- ❖ Study rapid size dynamics (~10sec res.)
- ❖ Interpretation of growth mechanisms
- ❖ Synthesis /formulation development

3. Larger scale process monitoring


Process development of drug-conjugated particles

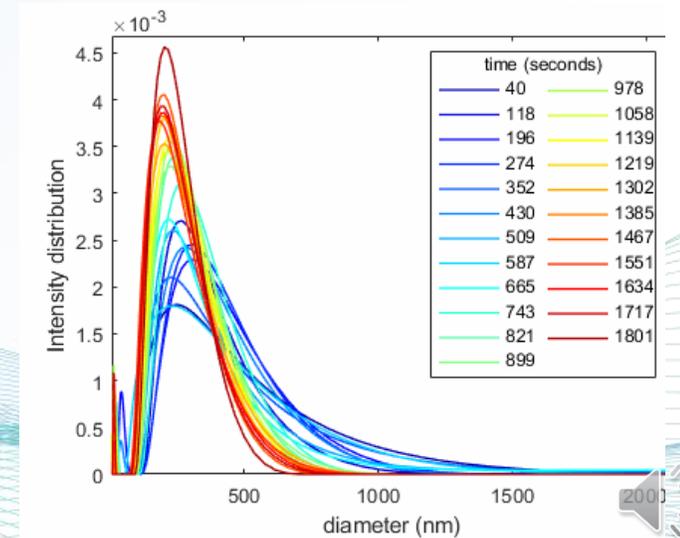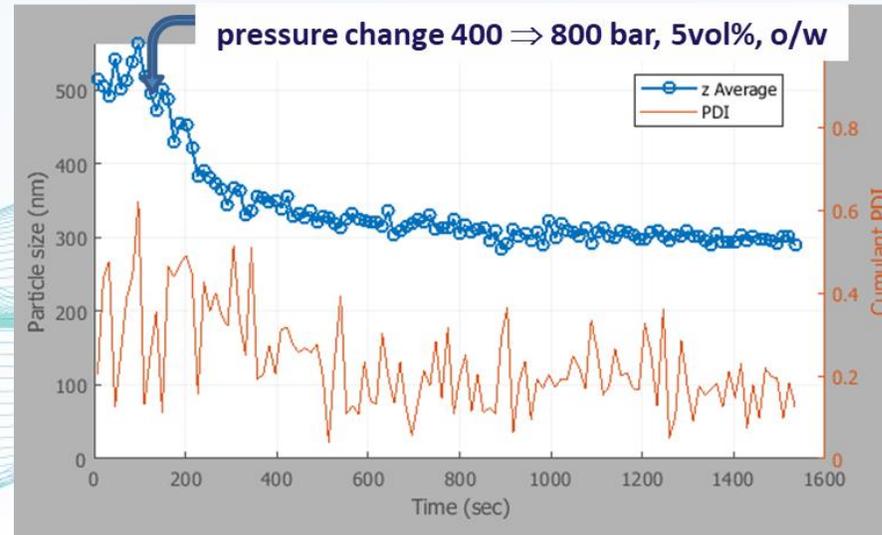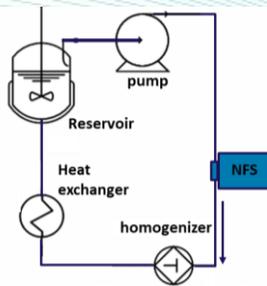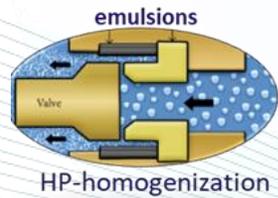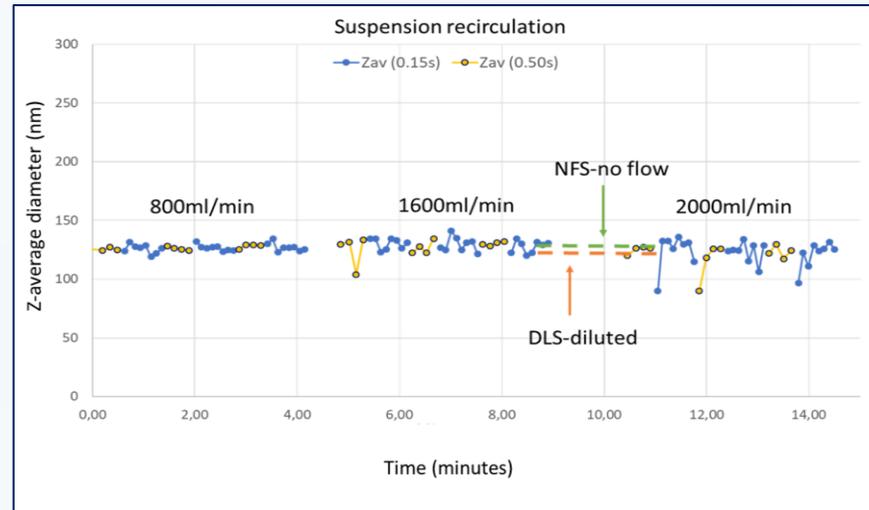

3. Application considerations

Flow requirements:

1. Non-turbulent, fully developed Poiseuille flow
2. Maximum ratio shear rate/diffusion rate (maximum 'Peclet' number)
 - both (1) and (2) are checked during measurement
 - upscaling via suitable flow cell / process connection

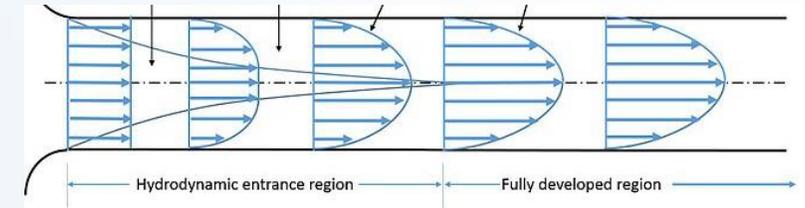

Concentration:

- High concentration => modifies diffusion & Stokes-Einstein relation
- Sizing possible (i) qualitatively or (ii) via calibration/knowledge of interactions
- We develop models (incl. hydrodynamics/structure) for quantitative sizing in this regime.

	Low Turbidity	High Turbidity
Low Concentration	✓	✓
High concentration	Calibration, modeling	Calibration, modeling

4. Summary

- **Spatially Resolved DLS: a significant step forward in nanosuspension characterization for Process Analytics and R&D**
 - **Rapid measurements of highly turbid and/or flowing samples for a broad range of sample /process environments.**
- => Process monitoring and/or Process control**



We are keen to explore new applications and further broaden the possibilities of the NanoFlowSizer.

Feel free to contact us to discuss your applications, request an 'NFS questionnaire' or our publication:

info@nanoflowsizer.com

<https://www.inprocess-lsp.com/contact/>

European Journal of Pharmaceutical Sciences 133 (2019) 205–213



Contents lists available at ScienceDirect

European Journal of Pharmaceutical Sciences

journal homepage: www.elsevier.com/locate/ejps

New unique PAT method and instrument for real-time inline size characterization of concentrated, flowing nanosuspensions

R. Besseling*, M. Damen, J. Wijgengangs, M. Hermes, G. Wynia, A. Gerich

InProcess-LSP, Kloosterstraat 9, 5349, AB, Oss, the Netherlands

Thank you for your interest !